

TECHNICAL NOTE

MesenCult™-ACF Plus Medium (Catalog #05445/05448)

Date of Notification: September 25, 2023

IMPORTANT: Please update your customers of this change using the following formal communication

I am writing to inform you that, as part of continual improvements to our products, we are **updating the recommended storage conditions for complete MesenCult™-ACF Plus Medium** (Catalog #05445/05448). You are being notified because you have ordered MesenCult™-ACF Plus Medium within the last 6 months.

The **new recommendation is as follows:** If not used immediately, complete MesenCult™-ACF Plus Medium (MesenCult™-ACF Plus Medium + MesenCult™-ACF Plus 500X Supplement + L-Glutamine) should be stored at 2 - 8°C for up to 1 week to ensure optimal performance.

We are working actively to update our product documentation to reflect this updated guidance. If you require more information about this update, please contact Product and Scientific Support at techsupport@stemcell.com.

If you are not the user of this kit, please forward this Product Notification to the appropriate person.