


# Navigator表示の変更

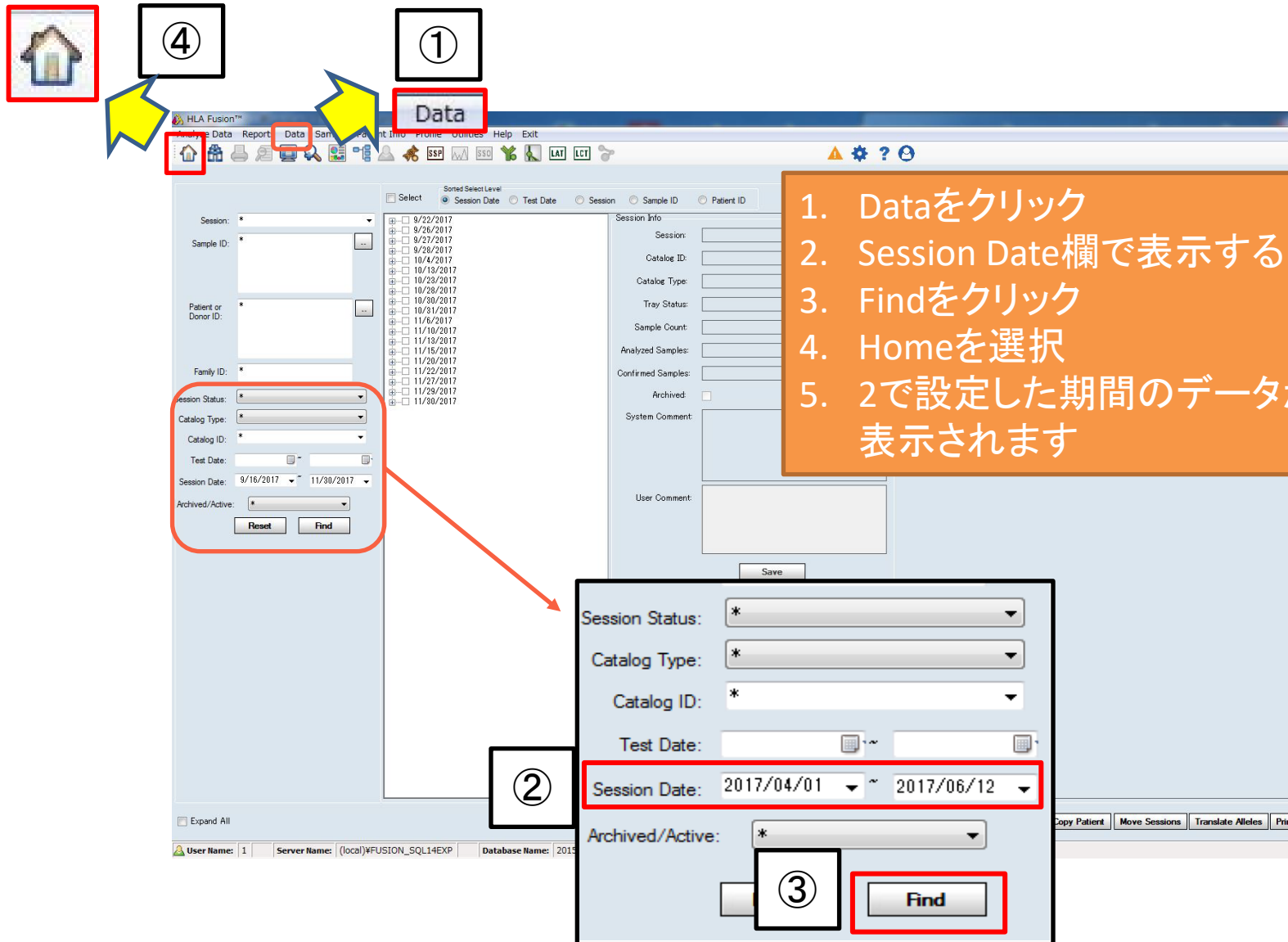

The screenshot shows the HLA Fusion software interface. The 'Data' menu item is highlighted with a red box and labeled '1'. The 'Home' icon in the toolbar is highlighted with a red box and labeled '4'. The 'Session Date' filter is highlighted with a red box and labeled '2', showing a date range from 2017/04/01 to 2017/06/12. The 'Find' button is highlighted with a red box and labeled '3'. A yellow arrow points from the 'Data' menu to the 'Home' icon. A red arrow points from the 'Session Date' filter to a magnified view of the filter controls.

1. Dataをクリック
2. Session Date欄で表示する日にちを設定
3. Findをクリック
4. Homeを選択
5. 2で設定した期間のデータがNavigatorに表示されます