


Navigator表示の変更

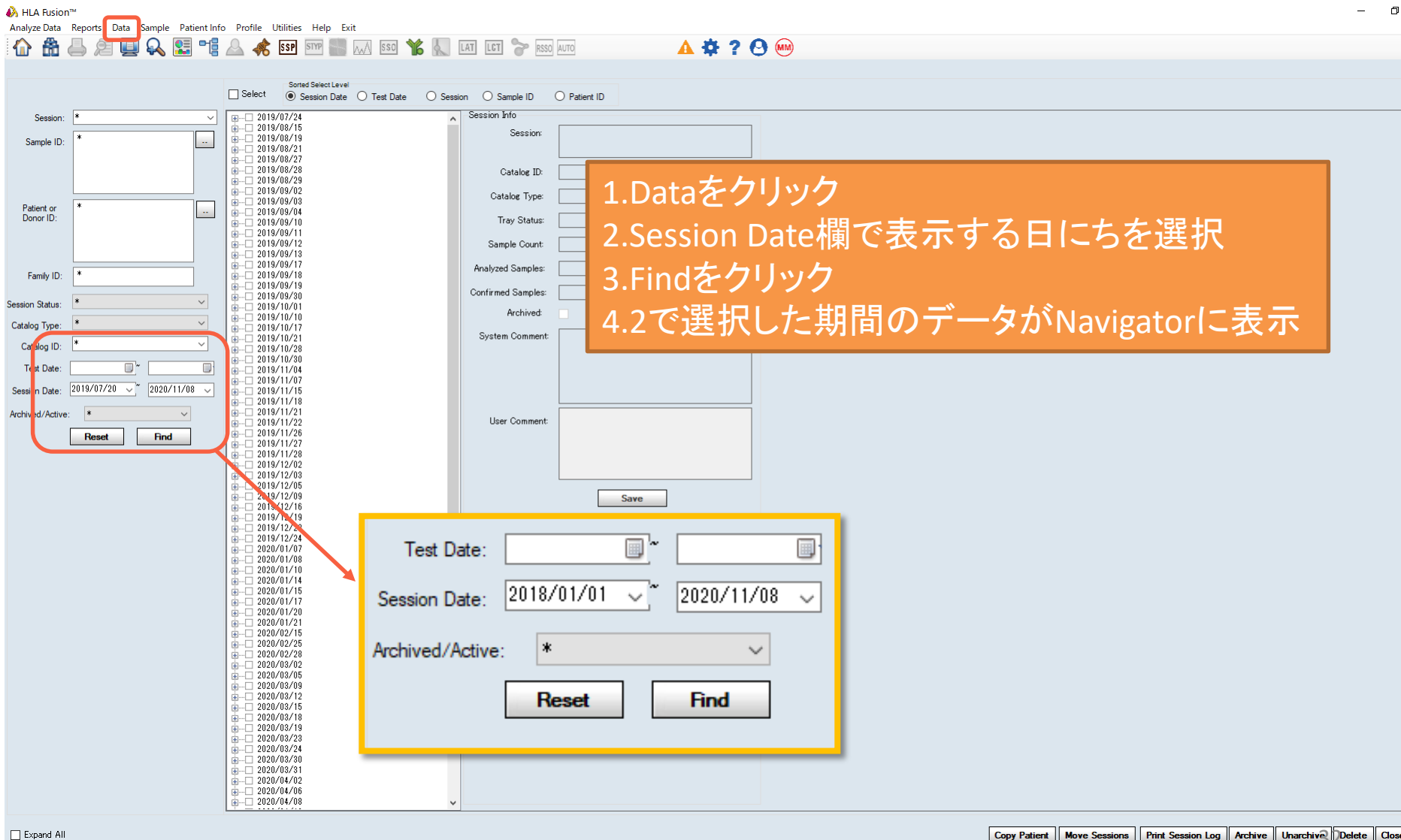

HLA Fusion™
Analyze Data Reports **Data** Sample Patient Info Profile Utilities Help Exit

Sorted Select Level
 Session Date Test Date Session Sample ID Patient ID

Session: *
Sample ID: *
Patient or Donor ID: *
Family ID: *
Session Status: *
Catalog Type: *
Catalog ID: *
Test Date: ~
Session Date: 2019/07/20 ~ 2020/11/08
Archived/Active: *

2019/07/24
2019/08/15
2019/08/19
2019/08/21
2019/08/27
2019/08/28
2019/08/29
2019/09/02
2019/09/03
2019/09/04
2019/09/10
2019/09/11
2019/09/12
2019/09/13
2019/09/17
2019/09/18
2019/09/19
2019/09/30
2019/10/01
2019/10/10
2019/10/17
2019/10/21
2019/10/28
2019/10/30
2019/11/04
2019/11/07
2019/11/15
2019/11/18
2019/11/21
2019/11/22
2019/11/26
2019/11/27
2019/11/28
2019/12/02
2019/12/03
2019/12/05
2019/12/09
2019/12/16
2019/12/19
2019/12/20
2019/12/24
2020/01/07
2020/01/08
2020/01/10
2020/01/14
2020/01/15
2020/01/17
2020/01/20
2020/01/21
2020/01/21
2020/02/15
2020/02/25
2020/02/28
2020/03/02
2020/03/05
2020/03/09
2020/03/12
2020/03/15
2020/03/18
2020/03/19
2020/03/23
2020/03/24
2020/03/30
2020/03/31
2020/04/02
2020/04/06
2020/04/08

Session Info
Session:
Catalog ID:
Catalog Type:
Tray Status:
Sample Count:
Analyzed Samples:
Confirmed Samples:
Archived:
System Comment:
User Comment:
Save

1. Dataをクリック
2. Session Date欄で表示する日にちを選択
3. Findをクリック
4. 2で選択した期間のデータがNavigatorに表示

Test Date: ~
Session Date: 2018/01/01 ~ 2020/11/08
Archived/Active: *
Reset Find

Expand All
Copy Patient Move Sessions Print Session Log Archive Unarchive Delete Close