



## CAUTION

The "EasyPlate" EasySep™ Magnet generates a strong magnetic field. Keep away from pacemakers, magnets, computer disks, watches, and other objects that respond to magnetic fields.

## THE "EASYPLATE" EASYSEP™ MAGNET IS ONLY FOR USE WITH CERTAIN EASYSEP™ HUMAN OR MOUSE ENRICHMENT KITS.

If your cell type or sample is not listed on [www.stemcell.com](http://www.stemcell.com), please contact us at [techsupport@stemcell.com](mailto:techsupport@stemcell.com).

## PRODUCT DESCRIPTION

A sealed anodized aluminum body housing rare earth magnetic elements.

## RECOMMENDED 96 WELL PLATE

The "EasyPlate" EasySep™ Magnet is designed to hold a 96 well plate (such as Costar, Catalog #3788 or BD Biosciences, Catalog #351177). Round bottom, non-tissue culture-treated plates work best.

If using a different type of non-tissue culture treated 96 well plate, ensure that it properly fits on the "EasyPlate" EasySep™ Magnet before use. Some 96 well plates may not sit flat on the magnet, which could affect the success of the separation.

## CARE OF MAGNET

After completing a cell separation, wipe the magnet clean with distilled water and then with alcohol to dry the magnet. Do not clean with bleach.

CATALOG #18102

Components:

- “EasyPlate” EasySep™ Magnet

 EasySep™

MAGNETS




WWW.STEMCELL.COM

FOR RESEARCH USE ONLY. NOT FOR THERAPEUTIC OR DIAGNOSTIC USE

AUGUST 2011

IN NORTH AMERICA TOLL-FREE T. 1 800 667 0322 TOLL-FREE F. 1 800 567 2899 T. +1 604 877 0713 F. +1 604 877 0704 E. INFO@STEMCELL.COM  
IN EUROPE TOLL-FREE T. 00 800 7836 2355 TOLL-FREE F. 00 800 7836 2300 E. INFO.EU@STEMCELL.COM  
IN AUSTRALIA TOLL-FREE T. 1 800 060 350 F. +61 (03) 9338 4320 E. INFO.AUS@STEMCELL.COM  
IN SINGAPORE T. +(65) 6776 7754 F. +(65) 6776 7114 E. INFO.SG@STEMCELL.COM

#28785  
V1.1.0

 Printed on recycled paper.